

# EYEWASH FLUSH LOG

**Turn on your emergency eyewash *at least* once a month for a minimum of 3 minutes and document in the log**

**Purpose of testing:** Confirm the eyewash works correctly, flush sediment from the supply line, and reduce microbial growth from standing water

✓ **PASS** criteria (after 3 minutes of operation)

- Water flow alone opens the covers without user assistance
- Water is clear and not too hot or cold
- Water reaches both eyes at the same time
- Streams of water maintain controlled flow and do not hurt the eyes



**✗ FAIL** if any pass criteria is not met or you have other concerns about the station functionality  
If FAIL: Submit a **work order** or email your **building manager**

- Include the room number
- Describe the eyewash location  
(*Example: across from bench #3.*)
- Describe what is not working

### Facilities Service Request:

BPSB, ERC/WERC, GCIS, Jones,  
Searle, Kent, Kersten (KPTC),  
Hinds, Hull Court (Zoology,  
Erman, Anatomy, Culver),  
Michelson (MCP)



**PLEASE EMAIL RESEARCHSAFETY@UCHICAGO.EDU FOR ANY ADDITIONAL CONCERNS**

[illegible][illegible][illegible][illegible]